

# Juno™ System Software

## v3.7.1

### v3.7.1 Enhancements

None.

### v3.7.1 Bug Fixes

**IFC orientation check.** Orientation for IFCs (integrated fluidic circuits) is now determined by the location of the barcode label. If the barcode is not detected, the user is prompted to check the orientation of the IFC.

## v3.5.1

### v3.5.1 Enhancement

**LP—48.48 IFC support.** New scripts to run the LP—48.48 IFC.

### v3.5.1 Bug Fixes

- **Leap year.** Scripts recognize 29 February as a valid date.
- **Daylight saving time.** Instrument clock adjusts for daylight saving time.

## v3.3.4

### v3.3.4 Enhancement

**Juno LP—192.24 IFC support.** New scripts to run the Juno LP—192.24 IFC.

### v3.3.4 Bug Fixes

None.

## v3.1.4

### v3.1.4 Enhancements

- **System power down.** This new operation on Juno permits safe shutdown of the operating system before powering down the instrument. While the system is running, press the front power and standby indicator for two seconds to initiate the shutdown.
- **Barcode recognition.** Improved to recognize all supported IFCs on the instrument.
- **System cleaning script.** Updated with a running time of 13 minutes.

### v3.1.4 Bug Fixes

None.

#### **For technical support visit [fluidigm.com/support](https://fluidigm.com/support).**

North America +1 650 266 6100 | Toll-free (US/CAN): 866 358 4354 | [techsupport@fluidigm.com](mailto:techsupport@fluidigm.com) Latin America +1 650 266 6100 | [techsupportlatam@fluidigm.com](mailto:techsupportlatam@fluidigm.com)  
Europe/Middle East/Africa/Russia +44 1223 859941 | [techsupporteurope@fluidigm.com](mailto:techsupporteurope@fluidigm.com) China (excluding Hong Kong) +86 21 3255 8368 | [techsupportchina@fluidigm.com](mailto:techsupportchina@fluidigm.com)  
Japan +81 3 3662 2150 | [techsupportjapan@fluidigm.com](mailto:techsupportjapan@fluidigm.com) All other Asian countries/India/Australia +1 650 266 6100 | [techsupportasia@fluidigm.com](mailto:techsupportasia@fluidigm.com)

---

#### **For Research Use Only. Not for use in diagnostic procedures.**

Information in this publication is subject to change without notice. **Patent and license information:** [fluidigm.com/legalnotices](https://fluidigm.com/legalnotices). Fluidigm, the Fluidigm logo, and Juno are trademarks or registered trademarks of Fluidigm Corporation in the United States and/or other countries. All other trademarks are the sole property of their respective owners. © 2017 Fluidigm Corporation. All rights reserved. 06/2017